

SoftGrip™ Pipette Bulletin

The New Hamilton SoftGrip™ Pipette

The new SoftGrip boasts an enlarged ejector button, dual needle volume indicator, slimmer ejector sleeve, redesigned plunger button, and a more stain resistant grip. These modifications have enhanced the functionality of the SoftGrip while maintaining the principals of quality, ergonomics, and functionality that made the SoftGrip popular.



SoftGrip™ Features and Benefits

- **Best overall accuracy and precision**
- **Fully Autoclavable**
- **Made in the USA**
- **Ergonomic by design**

Soft Contoured Handle

The SoftGrip handle reduces strain on the user resulting from continuous and repetitive operation.

Curved Hilt

The hilt allows the pipette to hang freely from the hand further reducing the hand strain associated with repetitive pipetting.

Low Plunger and Ejector Force

A soft plunger spring and enlarged ejector button minimizes the force required for operation. Decreased forces will result in less fatigue and repetitive stress injuries.

Color Coded Volumes

Unique color coding and labeling clearly indicate the volume of the pipette as well as the acceptable working range for adjustable volume pipettes.

Isolated Volume Adjustment

The unique location of the volume adjustment ring reduces accidental volume changes.

Improved Volume Indication

The dual needle indicator improves volume adjustment accuracy by removing parallax errors.

Universal Tips


The SoftGrip's universal tip nozzle ensures a proper fit with universal pipette tips. The quality of the tip used will affect the pipettes overall accuracy and precision. High quality Hamilton AdvanTips™ are recommended for all applications.

Slim Ejection Sleeve

The ejector sleeve is made from a high strength, chemically resistant polymer. The slim design allows the pipette to reach to the bottom of narrow tubes without the need for extended length tips.

Simple Calibration

A Universal Calibration Key reduces the need for costly calibration services. Calibration requires only minutes to perform and eliminates the downtime and cost of alternative calibration methods.

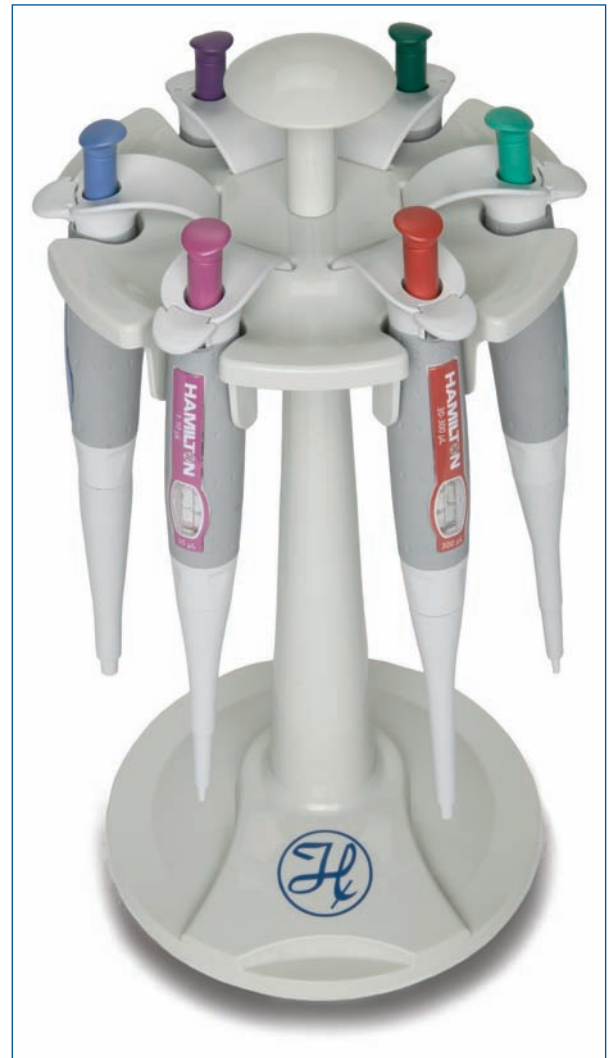

HAMILTON
THE MEASURE OF EXCELLENCE™

Hamilton Company
4970 Energy Way
Reno, Nevada 89520 USA
Toll Free: 800 648-5950
Telephone: +1-775-858-3000
Fax: +1-775-856-7259
e-mail: sales@hamiltoncompany.com

Hamilton Bonaduz AG
Via Crusch 8
CH-7402 Bonaduz/Switzerland
Toll-Free: 00800-660-660-60
Telephone: +41-(0)81-660-60-60
Fax: +41-(0)81-660-60-60
e-mail: marketing@hamilton.ch

www.hamiltoncompany.com